

## How to Use the MS Syringe Filters

### Warning

- Do not use the MS filter unit with male luer slip outlet for direct patient care applications; they are designed for laboratory use only.
- Do not use with syringes smaller than 10 cc because pressures in excess of the maximum pressure rating may be reached, potentially causing damage to the filter unit and/or personal injury.

### Cautions

- Do not use the MS syringe filters at temperatures above 45°C (113 °F) to ensure proper filtration.
- Do not use the MS unit to filter a protein containing solution unless binding of the sample to the filter is evaluated.
- Do not reuse the MS 2 filter unit with male luer slip outlet.

### Directions for the MS Filter Unit with Male Luer Slip Outlet


1



•

Fill the syringe with the solution you want to filter.

2

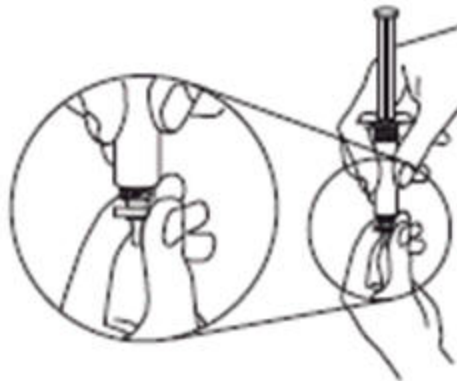

•


• Attach the syringe to the MS filter unit with male luer slip outlet.

3



- 
- Hold the syringe with filter pointing up and "top off" by pushing a few drops through the filter.

4

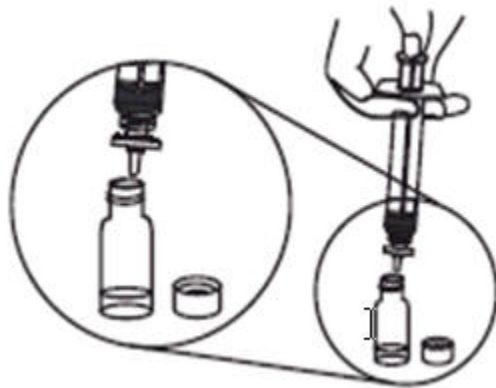

- 
- Push the syringe plunger to deliver a filtered solution.